



LUMIGEN, INC.
A Beckman Coulter Company

22900 W. Eight Mile Rd., Southfield, MI 48033
Tel (248) 351-5600 Fax (248) 351-0518

Lumigen[®]
Lumi-Phos[®] 530 & Lumi-Phos[®] Plus
Suggested Product Handling Instructions

Storage:

Store bottled product at 2-8° C.
Protect from light contamination.

Do not freeze.

Use:

*Use in subdued light. Indirect incandescent lighting is preferred.
Exposure to direct light will cause elevated background.

- 1) Allow solution to come to room temperature (Approximately 1 hour for 100 mL of product).
- 2) Gently invert (4-5 times) solution in its packaged containers to assure homogeneity prior to dispensing.
- 3) Dispense the amount needed of solution into a new container. Containers should be opaque or covered with aluminum foil to protect from direct light (daylight or artificial).
- 4) **Do not mix again before use.**
- 5) Aliquot the working solution into assay wells or gel, while protecting from light contamination.
- 6) Store unused product in original container at 2-8° C for up to 2 years.

Repackaging:

Repackaging of Lumigen products is discouraged as reliability can be compromised by contamination.

Bottling is available from Lumigen to suit the volume demands for your specific work process.

If you choose to repackage, new opaque HDPE or PP plastic containers are required. The high temperatures in the container molding process destroy potential contaminants. **Reusing and washing of containers can lead to contamination and subsequent high background.**